
Dragon 200 Lab Burner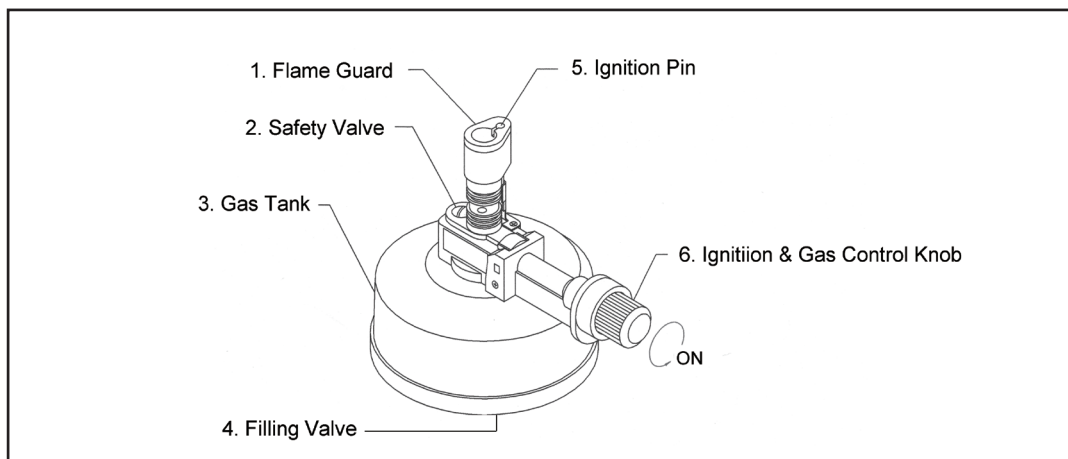

ATTENTION: This unit does not contain any butane gas when purchased. Please follow below to make it operational.

1. FUEL FILLING AND REFILLING

- Make sure the gas burner is switched off before filling.
- Reverse the unit upside down.
- Insert the container's nozzle vertically in a downward motion into the FILLING VALVE ④ on the bottom of the unit.
- When the fuel begins to overflow, remove the nozzle from the filling valve to stop the filling operation. It means the unit is full of gas.
- Allow a few minutes for the fuel to stabilize in the fuel tank.

2. IGNITION

- Turn the IGNITION BUTTON ⑥ to the left (counterclockwise) about 90 degree.
- Press the IGNITION BUTTON ⑥ to light the flame.

3. SHUT-OFF

Turn the IGNITION BUTTON ⑥ to right (clockwise) to decrease until the end to shut off the flame.

4. FLAME LENGTH ADJUSTMENT

After the ignition, turn the GAS CONTROL KNOB ⑥ to the left (counterclockwise) to increase the flame; to the right (clockwise) to decrease the flame and shut if off.

4. WARNING

- Do not touch FLAME GUARD ① when hot.
- Keep out of reach of children.
- Do not drop. Do not operate or store in temperature exceeding 50°C.
- Do not try to disassemble the unit.

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