

## Rocker 300 - MF 31 Vacuum Filtration System

Vacuum filtration is widely applied in laboratory for various tests such as microorganism test in food, pharmacy, beverage and drinking water in water treatment field.

Rocker 300-MF31 is widely used in suspended solid test and general filtration.

### Unpacking

When unpacking, make sure the below items are included.

1. Vacuum Pump Rocker 300... .....1 ea.
2. MF31, PES Filtration Set... ..1 ea.
3. Silicone tube (1m)... ..1 ea.



### Assembling the system

1. Cut the silicon tube into suitable length.
2. Connect the tube to the air inlet of the pump and the other to the nozzle of the glass set.

### Ordering information

1. 167311-11 Vacuum filtration system Rocker 300 - MF 31 · AC 110V/60Hz
2. 167311-22 Vacuum filtration system Rocker 300 - MF 31 · AC 220V/50Hz
3. 167300-11 Oil free vacuum pump Rocker 300, AC 110V/60Hz
4. 167300-22 Oil free vacuum pump Rocker 300, AC 220V/50Hz
5. 200320-31 MF31, PES Filtration Set
6. 200300-01 MF3, Magnetic Filter Holder 300ml with lid kit
7. 167100-15 Receiver Flask 1000ml with drain and silicone fixing sucker
8. 167200-38 Silicon tube (1m)
9. 167100-20 PALL, Type A/ E Glass Fiber Discs, 47mm, 1 um, 100/PK
10. 167100-52 PALL, GN-6 Metrical Membrane Disc Filter, grid, 47mm, 0.45 um, 200/PK