


VF10 Glass Filtration Set

Unpacking

Carefully check if the shipping carton is damaged before unpacking. Please check all supplied items according to the standard package list showed as the right photo. If you have any questions, please contact your local dealer and provide with serial number & shipping carton for help.

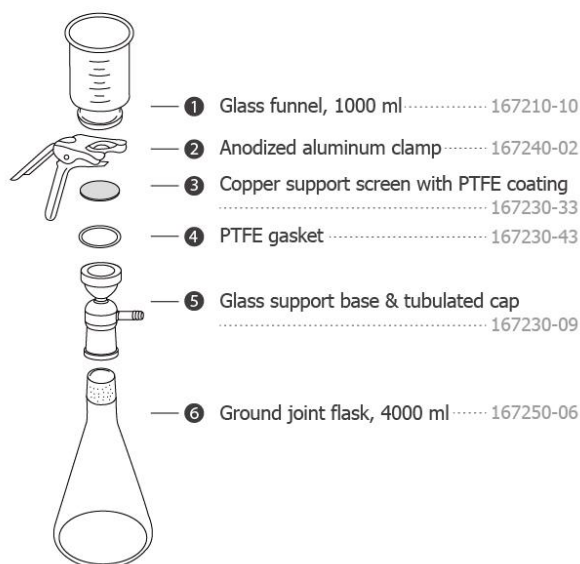

Usage Guide

1. Put the glass support base into the flask.
2. Put the PTFE gasket and the copper support screen into the support base then put 47mm membrane filter on the screen.
3. Put the glass funnel on the glass support base, and clamp them together.

***** Please confirm all steps above are completed before doing filtration test.**

Ordering Information

167200-10:	VF10 Glass filtration set
167210-10:	1000ml Glass funnel (90mm)
167240-02:	Anodized aluminum clamp
167230-33:	90mm Copper support screen with PTFE coating
167230-43:	90mm PTFE gasket
167230-09:	90mm Glass support base & tubulated cap
167250-06:	4000ml Ground joint flask