


## VF11 Stainless Steel Filtration Set

### Unpacking

Carefully check if the shipping carton is damaged before unpacking. Please check all supplied items according to the standard package list showed as the right photo. If you have any questions, please contact your local dealer and provide with serial number & shipping carton for help.

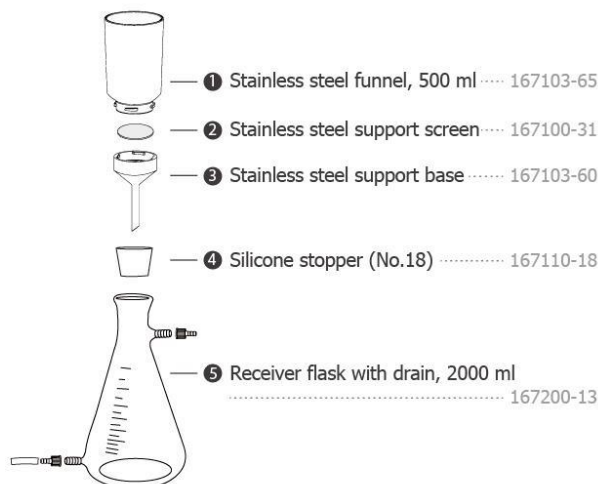

### Usage Guide

1. Put the fixing sucker on a flat table.
2. Press down the fixing sucker from the middle to squeeze the air outward, and make sure the "holding base" is connected to the table tightly.  
( It is easier to remove the fixing sucker from table by taking the "Handle" )
3. Put the glass flask on the "Holding base" and let the "Holding side" cover the glass flask.
4. Put the SS support base and silicone stopper together, and connect to the flask.
5. Screw the screwthread connectors onto the hose barb on flask, then connect a matched tube to drain hose barb.
6. Put the SS support screen into the support base then put 47mm membrane filter on the SS support screen.
7. Put the SS funnel on the SS support base then spin to lock them together.

**\*\*\* Please confirm all steps above are completed before doing filtration test.**

### Ordering Information

167200-11:	VF11 SS filtration set
167103-65:	500ml SS funnel (47mm)
167100-31:	47mm SS support screen
167103-60:	47mm SS support base
167110-18:	No. 18 Silicone stopper
167200-13:	2000ml Receiver flask with drain