

## VF15 Glass Filtration Set

### Unpacking

Carefully check if the shipping carton is damaged before unpacking. Please check all supplied items according to the standard package list showed as the right photo. If you have any questions, please contact your local dealer and provide with serial number & shipping carton for help.

### Important Notice

- The product is suitable for GL45 storage bottle and fi
- Made of PP material filtration adaptor resistant to high temperature up to 130°C and autoclavable.
- Do not sterilize the product with flame.

### Usage Guide

1. Please ensure that 『 PTFE Gasket 』 has been installed properly in the filtration adaptor.
2. Screw the adaptor onto storage bottle tightly.
3. Put the glass support base and silicone stopper together, and connect to storage bottle by placing No. 8 stopper on the upper receiving port of adaptor.
4. Ensure the PTFE gasket and support screen has been installed into the support base then put membrane filter on the SS support screen.
5. Put the glass funnel on the glass support base then clamp them together.
6. Connect a matched tube between its hose barb and vacuum source to start filtration.

### Ordering Information

167200-15:	VF15 Glass filtration set
167210-10:	1000ml Glass funnel (90mm)
167240-02:	Anodized aluminum clamp
167230-33:	90mm PTFE coating membrane support
167230-43:	90mm PTFE gasket
167230-10:	90mm Glass support base
197000-65:	Filtration adaptor
197000-55:	5000ml Storage bottle (Duran with GL45)

