


## VF3 Glass Filtration Set

### Unpacking

Carefully check if the shipping carton is damaged before unpacking. Please check all supplied items according to the standard package list showed as the right photo. If you have any questions, please contact your local dealer and provide with serial number & shipping carton for help.

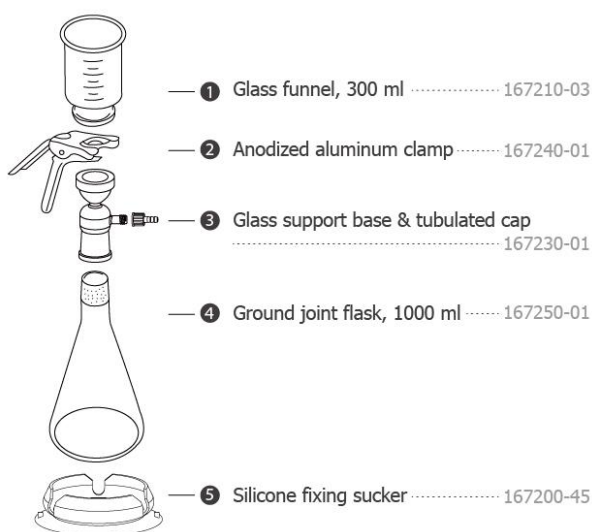

### Usage Guide

1. Put the fixing sucker on a flat table.
  2. Press down the fixing sucker from the middle to squeeze the air outward, and make sure the "holding base" is connected to the table tightly.  
( It is easier to remove the fixing sucker from table by taking the "Handle" )
  3. Put the flask on the "Holding base" and let the "Holding side" cover the glass flask.
  4. Screw the screwthread connector onto the hose barb then connect the glass support base to the flask.
  5. Put a 47mm membrane on the glass support base.
  6. Put the glass funnel on the glass support base then clamp them together.
- \*\*\* Please confirm all steps above are completed before doing filtration test.**

### Ordering Information

|            |   |
|------------|---|
| 167200-03: | VF3 Glass filtration set                  |
| 167210-03: | 300ml Glass funnel (47mm)                 |
| 167240-01: | Anodized aluminum clamp                   |
| 167230-01: | 47mm Glass support base & tubulated cap   |
| 167250-01: | 1000ml Ground joint flask                 |
| 167200-45: | Silicone fixing sucker (for 1000ml flask) |