


Rocker's multilayer filtration adaptor can be used with 2 kinds of membrane at the same time to improve the efficiency and reduce the repetitive work. The multilayer support base is available to connect with VF 3, VF 6, VF 7, and VF 12., the 47 mm glass filtration apparatus.

Carefully check if the shipping carton is damaged before unpacking. Please check the items according to package list. If you have any question, please preserve serial no. & shipping carton and contact your dealer for quick solution.

Item	Assembly
1	47 mm SS Support Screen
2	47 mm PTFE Gasket
3	47 mm Multilayer Support Base

Note. An additional clamp is required when set up with a multilayer support base.

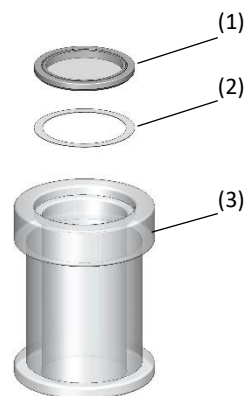

Assembling

Assembling (Please follow the diagram to install the multilayer support base)

1. Assemble the multilayer filtration adaptor according to above diagram.
2. Install the assembled adaptor between funnel and support base.
3. Place the desired membrane on each support base and fixed it with aluminum clamp.
4. Pour the liquid, turn on the pump and start a filtration.



Maintenance

1. Please make sure the pore size of upper layer membrane is bigger than down layer to ensure the fluently filtration.
2. To prevent air leakage, please make sure clamp all components right and tight.

Ordering Information

167230-05	Multilayer Filtration Adaptor	167230-44	Multilayer Support Base
167230-31	47 mm SS Support Screen	167230-41	47 mm PTFE Gasket
167240-01	Aluminum clamp	167200-07	VF 7, Glass Filtration Set